

Method 508 mod RE 2/11/02 Page 78
re-exts Ref HGMS-16965

Lab #	Description	Sample vol/wt	Drip rate ml/min	Surr add	Sol add	IS add	Final vol.	T.V. pos.	pH	Count
LB		500 mL		1 mL			1 mL		7	
Ref					0 mL					
Ref										
31662								9A		
31663								2B		
00006		460 mL						3A		
30643	has ref	500 mL						2A		
30645	2							1B		
30647	2							1A		
31460								3B		
Remarks 30645 and 30647 were brought to a pH of 7 by adding 1 1/2 mL of NaOH @ 10M to each										
drip rate must be between 5-15 ml/min for liq-liq extractors drip rate -										
Extraction type - <u>sep fun</u>		Solv. lot# <u>V41470</u>		NaCl lot# <u>V39618</u>		Na2SO4 lot# <u>V391400</u>				
STANDARD	INT STD LOG#									
CALIBRATION	<u>11201300</u>									
REF/SPIKE	<u>91759</u>									
MDL/MDRF										
Q. CHECK										
INT. STD.	<u>1561300</u>									
SURROGATE	<u>020102</u>									
LPC										
ENDRIN/DDT										

Reviewed by [Signature]
 Date 6/6/02

Analyst LP
 revision 4/12/01

Verified by [Signature]

Reviewed by [Signature]
 Date 7/15/02

Sequence Name: C:\HPCHEM\1\SEQUENCE\FEB1302.S

Comment: INST 209; 30Mx0.25umX0.50

Operator: 209 GALOS

Data Path: C:\HPCHEM\1\data\feb1302\

Pre-Seq Cmd:

Post-Seq Cmd:

Method Sections To Run

☒ Full Method☐ Reprocessing Only

On A Barcode Mismatch

☒ Inject Anyway☐ Don't Inject

Line	Type	Vial	DataFile	Method	Sample Name
1	Sample	48	940	GCMS0208	GC/MS SCAN 1/14
2	Sample	49	941	GCMS0208	GC/MS SCAN 1/14
3	Sample	50	942	GCMS0208	GC/MS SCAN 1/14
4	Sample	3	LB	GCMS0208	GC/MS SCAN 2/11
5	Sample	4	REF	GCMS0208	GC/MS SCAN 2/11
6	Sample	5	REFA	GCMS0208	GC/MS SCAN 2/11
7	Sample	6	31662	GCMS0208	GC/MS SCAN 2/11
8	Sample	7	31663	GCMS0208	GC/MS SCAN 2/11
9	Sample	8	6	GCMS0208	GC/MS SCAN 2/11
10	Sample	9	30643	GCMS0208	GC/MS SCAN 2/11
11	Sample	10	30645	GCMS0208	GC/MS SCAN 2/11
12	Sample	11	30647	GCMS0208	GC/MS SCAN 2/11
13	Sample	12	31460	GCMS0208	GC/MS SCAN 2/11
14	Sample	13	30644	GCMS0208	GC/MS SCAN
15	Sample	14	30648	GCMS0208	GC/MS SCAN
16	Sample	15	31133	GCMS0208	GC/MS SCAN

DFTPP

Data File : C:\HPCHEM\1\DATA\FEB1102\DFTPP3.D

Acq On : 13 Feb 2002 8:53 am

Sample :

Misc :

MS Integration Params: GOOFY.P

Method : C:\HPCHEM\1\METHODS\GCMS0208.M (RTE Integrator)


Title : 209 625/TCLP calibration table

Vial: 1

Operator: 209 GALOS

Inst : GC/MS 209

Multiplr: 1.00

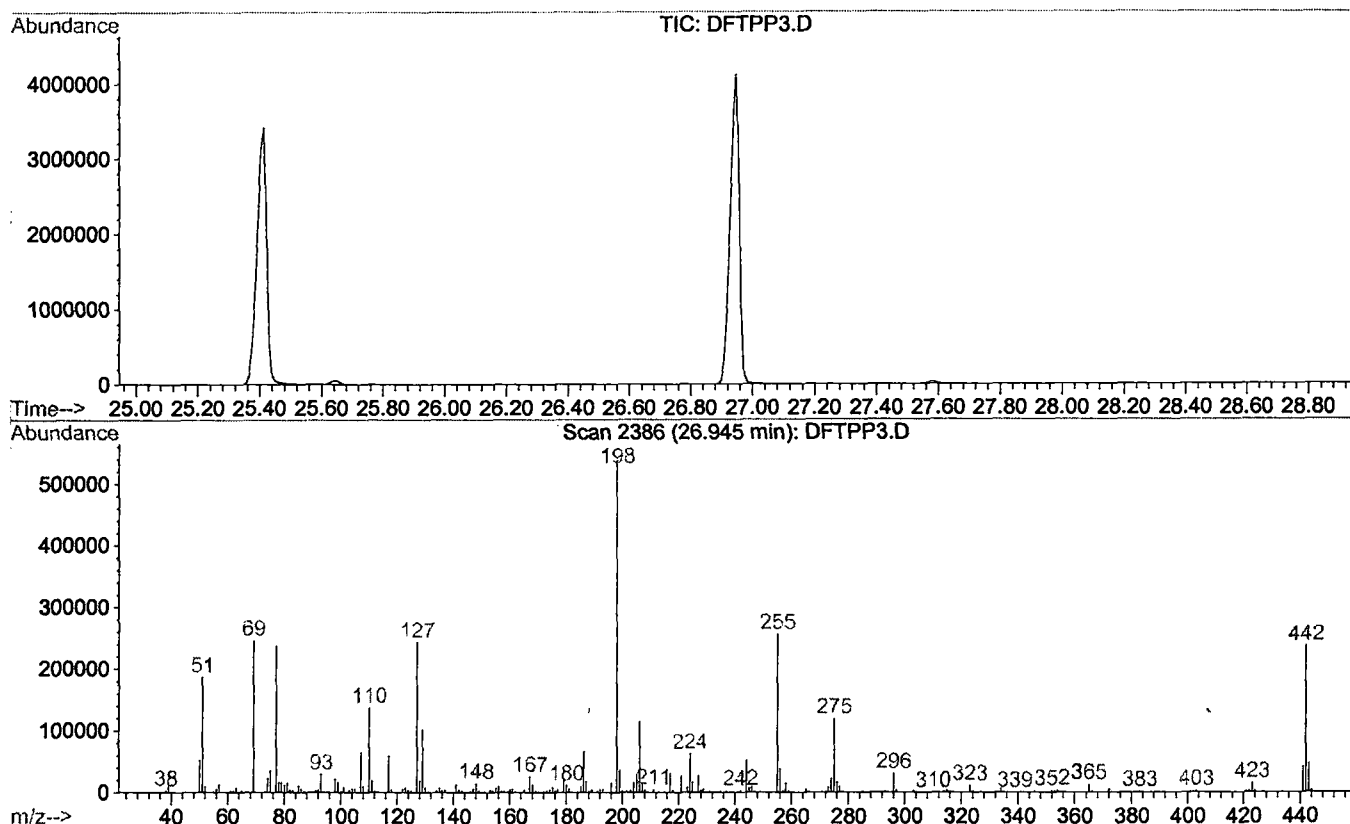

Spectrum Information: Scan 2386

Target Mass	Rel. to Mass	Lower Limit%	Upper Limit%	Rel. Abn%	Raw Abn	Result Pass/Fail
51	198	30	60	34.8	186944	PASS
68	69	0.00	2	0.0	0	PASS
69	198	0.00	100	45.8	246272	PASS
70	69	0.00	2	0.0	0	PASS
127	198	40	60	45.3	243264	PASS
197	198	0.00	1	0.0	0	PASS
198	198	100	100	100.0	537408	PASS
199	198	5	9	6.6	35704	PASS
275	198	10	30	22.1	118800	PASS
365	198	1	100	2.3	12594	PASS
441	443	0.01	100	90.5	41408	PASS
442	198	40	110	43.9	236096	PASS
443	442	17	23	19.4	45776	PASS