

Method

508 MOD.

Date Extracted

4/3/02

Page

42

Rob

\$GCMS - 17957

Lab #	Description	Sample vol/wt	Drip rate ml/min	Surr add	Spk add	IS add	Final vol.	T.V. pos.	pH	Count		
L.B.		500 ml		1 mL			1 mL	9A	7			
Ref 1					1 mL			9A				
Ref 2					1 mL			8A				
4945	Re-EXT.							9B				
7633								7B				
7634	F.B.							10A				
7635								10B				
7636								8B				
7637								8B				
7638	FB							7B				
7765								8A				
5286	RE-EXT. See 4/12	↓		↓			↓	7A	↓			
Remarks												
Reviewed by 4/5/6/02												
drip rate must be between 5-15 ml/min for liq-liq extractors												
drip rate -												
Extraction type - SEP. FUN. Solv. lot# X01E19 NaCl lot# V39614 Na2SO4 lot# V46623												
STANDARD	INT STD LOG#											
CALIBRATION												
REF/SPIKE	2193											
MDL/MDRF												
Q. CHECK												
INT. STD.												
SURROGATE												
LPC												
ENDRIN/DDT												

Analyst

A.D.

Revision 4/12/01

Verified by

Reviewed by

Date

Dr
4/18/02

Sequence Name: C:\MSDchem\1\sequence\020410.S

Comment: GCMS scan

Operator: Nefliu

Data Path: C:\MSDchem\1\DATA\020410\

Pre-Seq Cmd:

Post-Seq Cmd:

Method Sections To Run On A Barcode Mismatch

(X) Full Method (X) Inject Anyway

() Reprocessing Only () Don't Inject

Line	Type	Vial	DataFile	Method	Sample Name
------	------	------	----------	--------	-------------


1	Sample	1	DFTPP	508FAST	dftpp
2	Sample	1	DFTPP2	508FAST	dftpp
3	Sample	2	0410C3	508SCAN1	ccv
4	Sample	3	0403LB	508SCAN1	04/03 lab blk
5	Sample	4	0404LB	508SCAN1	04/04 lab blk
6	Sample	5	0403REF1	508SCAN1	04/03 ref
7	Sample	6	0404REF1	508SCAN1	04/04 ref
8	Sample	7	0403REF2	508SCAN1	04/03 ref
9	Sample	8	0404REF2	508SCAN1	04/04 ref
10	Sample	9	07634	508SCAN1	field blk
11	Sample	10	07638	508SCAN1	field blk
12	Sample	11	04945	508SCAN1	sample
13	Sample	12	05286	508SCAN1	sample
14	Sample	13	07633	508SCAN1	sample
15	Sample	14	07635	508SCAN1	sample
16	Sample	15	07636	508SCAN1	sample
17	Sample	16	07637	508SCAN1	sample
18	Sample	17	07765	508SCAN1	sample
19	Sample	18	08019	508SCAN1	field blk
20	Sample	19	08012	508SCAN1	sample
21	Sample	20	08013	508SCAN1	sample
22	Sample	21	08014	508SCAN1	sample
23	Sample	22	08015	508SCAN1	field blk
24	Sample	23	08016	508SCAN1	sample
25	Sample	24	08017	508SCAN1	sample
26	Sample	25	08018	508SCAN1	sample

↓ ignore stopped here.

Method 508 GC MS scan
Instrument 214

File Directory 020410

Storage Directory _____

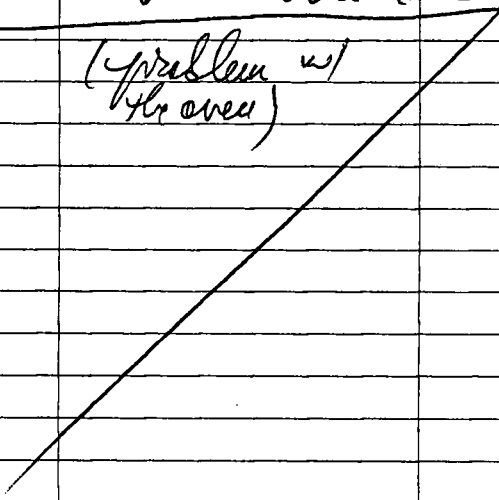Lab #	Filename	Description	Vial #		Inj.	Standards used - Lot #			Remarks	Initials
			front	rear		Calibration	Reference	Surrogate		
	4FTPP	dftpp	1		1					AN
	4FTPP2		1			2				
	041003	cen	2			240/2964				1
	0403LB	ens, bk	3							
	0404LB		4							
	0403Ref1	ref	5				293			
	0404Ref1		6							
	0403Ref2		7							
	0404Ref2		8							
	07634	sample	9							
	07638		10							
	04945		11							
	05286		12							
	07633		13							
	07635		14							
 (problem w/ the oven)										

Date 04/10/02


Extraction date 04/05/04/02

Page 3

DFTPP

Data File : C:\MSDCHEM\1\DATA\020410\DFTPP2.D
 Acq On : 10 Apr 2002 10:15 am
 Sample : dftpp
 Misc :
 MS Integration Params: EVENTS.E
 Method : C:\MSDCHEM\1\METHODS\SCANAPR0902.M (Chemstation Integrator)
 Title : Method 508 GC/MS Scan

Vial: 1
 Operator: Nefliu
 Inst : Instrume
 Multiplr: 1.00

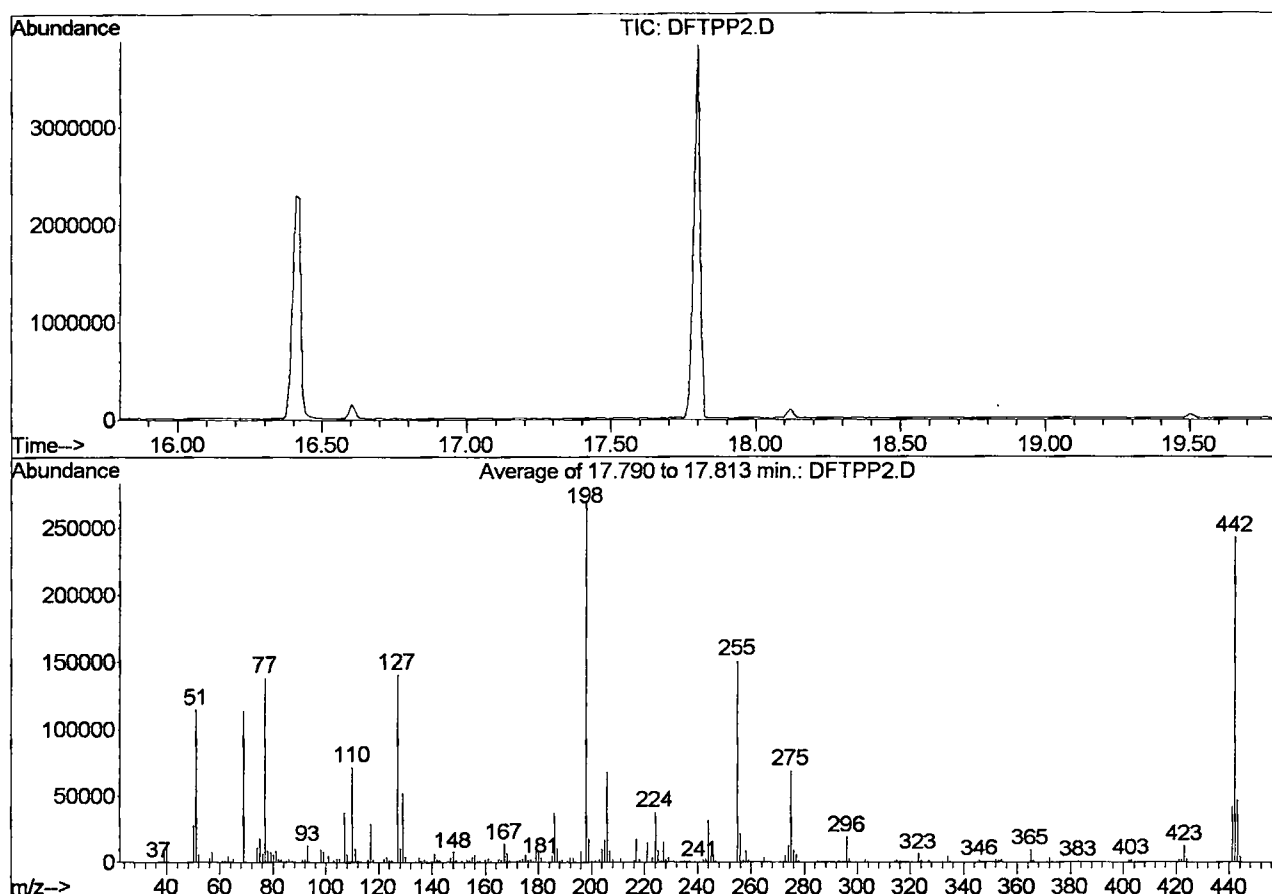

Spectrum Information: Average of 17.790 to 17.813 min.

Target Mass	Rel. to Mass	Lower Limit%	Upper Limit%	Rel. Abn%	Raw Abn	Result Pass/Fail
51	198	30	60	42.5	114808	PASS
68	69	0.00	2	0.0	0	PASS
69	198	0.00	100	42.3	114394	PASS
70	69	0.00	2	0.5	542	PASS
127	198	40	60	52.3	141344	PASS
197	198	0.00	1	0.0	0	PASS
198	198	100	100	100.0	270266	PASS
199	198	5	9	6.8	18284	PASS
275	198	10	30	25.6	69144	PASS
365	198	1	99	3.5	9531	PASS
441	443	1	100	89.8	42600	PASS
442	198	40	110	90.0	243298	PASS
443	442	17	23	19.5	47437	PASS